

DepleteX™ Single Cell RNA Boost Kit (Human) with compatible panels

KIT1018, KIT1024, KIT1025

Contents

Product overview.....	2	Protocol.....	8
Workflow.....	3	Step A: Ribonucleoprotein complex formation.....	8
Kit contents and storage.....	4	Step B: CRISPR digestion.....	10
Required materials and equipment provided by the user.....	6	Step C: Size Selection.....	12
Best Practices.....	6	Library validation.....	13
Revision log.....	7	Bioinformatics guidelines.....	14

Contact Us

If you have any questions, contact Technical Support at support@jumpcodegenomics.com

Find us at our website: jumpcodegenomics.com

Call us at 1.619.900.1701

Product overview

Traditionally, single cell data processing incorporates certain filtering and normalization steps prior to canonical clustering and downstream interpretation. Instead of filtering those reads in silico, DepleteX™ removes those reads in vitro, redistributing sequencing reads to biologically relevant transcripts.

Jumpcode depletion technology leverages CRISPR and a specifically designed guide RNA set to remove reads that are not incorporated in downstream analysis. The DepleteX Single Cell RNA Boost Kit gives you the ability to cut through the noise, with minimal impact on your workflow, and maximize your view of biology.

Content for depletion was designed by analyzing a cohort of publicly available 10x Genomics® Chromium single cell datasets from various sources. CRISPR complexes are combined with cDNA to deplete highly expressed protein-coding ribosomal and mitochondrial genes, non-variable genes that vary little in expression between different tissue types, and reads that do not align to the transcriptome. A list of the intervals targeted for depletion is provided on the Jumpcode Genomics website.

Depletion content is provided as two separate panels of 1) base content (ribo/mito) and non-transcriptomic intervals and 2) non-variable genes. Users have the flexibility to deplete or not deplete targeted non-variable genes.

DepleteX™ Single Cell RNA Boost Kit (Human)

Description	
Assay time	2 hours
Hands-on time	45 min
Input	Uses one of four Chromium cDNA aliquots per prep
Method	Single cell 3' gene expression libraries for 10x Genomics
10X Genomics Compatible Partner Program verified	10x Genomics Chromium Next GEM Single Cell 3' Reagent Kit v3.1

Jumpcode compatible products for use in this protocol:

Compatible Jumpcode products for this user manual	Depletion content
KIT1018 DepleteX™ Single Cell RNA Boost Kit (Human)	Human unaligned reads, ribosomal, mitochondrial, and non-variable genes
KIT1024 RNA Depletion Panel (Globin)	Human globin genes: HBA1, HBA2, HBB, HBD
KIT1025 RNA Depletion Panel (Insulin)	Human insulin gene: INS

Workflow

DepleteX™ Single Cell RNA Boost is a simple 3-step protocol easily integrated into 10x Genomics Chromium Next GEM Single Cell 3' gene expression protocol. This workflow can also be adapted to any single cell or spatial method for depleting undesired fragments from cDNA prior to sequencing.

The user follows the Chromium Next GEM Single Cell 3' protocol v3.1, without modification, until the end of **Step 3.4 – Post Ligation Cleanup – SPRIselect**. Immediately after Step 3.4, the user switches to the DepleteX™ protocol. After the depletion protocol is complete, the user resumes the Chromium Next GEM Single Cell 3' protocol v3.1 at **Step 3.5 – Sample Index PCR** and follows the protocol until completion of the libraries.

Protocol Overview:

- A. Perform library preparation up to and including Step 3.4 post-ligation cleanup
 - Step 1: GEM generation and barcoding
 - Step 2: Post GEM-RT cleanup and cDNA amplification
 - Step 3: Gene expression library construction
 - At the end of **Step 3.4**, elute the sample in 16 µL of Nuclease-Free Water (instead of 30.5 µL of Buffer EB).
- B. **DepleteX Single Cell RNA Boost Kit**
 - Ribonucleoprotein complex formation
 - CRISPR digestion
 - Size selection
- C. Continue with Chromium Next GEM Single Cell 3' protocol until completion of libraries
 - Step 3.5: Sample Index PCR
 - Step 3.6: Post Sample Index PCR Double Sided Size Selection - SPRIselect

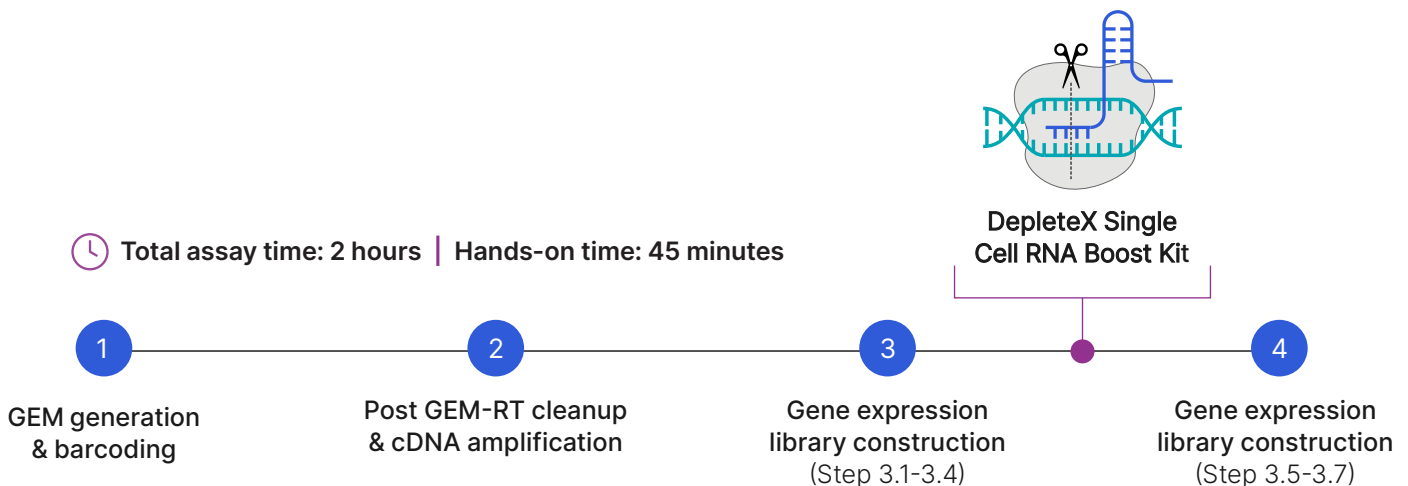


Figure 1: Schematic illustrating the position where DepleteX protocol fits into 10x Genomics Chromium Next GEM Single Cell 3' library preparation protocol.

Kit contents and storage

DepleteX™ Single Cell RNA Boost Kit (Human) contains enough material to deplete 24 Chromium single cell 3' gene expression libraries. The products listed below are sold separately. Kit contents and storage temperatures are indicated in the tables below.

Compatible Jumpcode products for use in this protocol:

- KIT1018 DepleteX Single Cell RNA Boost Kit (Human): ASY-1060, ASY-1061
- KIT1024 RNA Depletion Panel (Globin): ASY1064, ASY1071
- KIT1025 RNA Depletion Panel (Insulin): ASY1064, ASY1072

ASY-1060: DepleteX™ Single Cell RNA Boost Kit depletion reagents

24 samples

Stored at -20°C

Kit contents	Part number	Quantity
Cas9	REA1000	1
10x Cas9 Buffer	REA1001	1
RNase Inhibitor	REA1007	1
Nuclease-free Water*	REA1023	2

* **Note:** Store at 4°C after first use.

ASY-1061: DepleteX™ Single Cell RNA Boost Guide RNA

24 samples

Stored at -80°C

Kit contents	Part number	Quantity
Guide RNA Single Cell RNA - Boost	REA1025-003	1
Guide RNA Single Cell RNA - NVG	REA1120-001	1

ASY-1064: Depletion reagents

8 samples per box

Stored at -20°C

Kit contents	Part number	Quantity
Cas9	REA1039	1
10x Cas9 Buffer	REA1040	1
RNase Inhibitor	REA1041	1
Nuclease-Free Water	REA1042	1
Tris Buffer	REA1043	1
P5 Primer	REA1044	1
P7 Primer	REA1045	1
2X PCR Mix	REA1046	1

ASY-1071 Guide RNA for Globin

24 samples per tube

Kit contents	Part number	Quantity	Storage temperature
Guide RNA for Globin	REA1053	1	-80°C

ASY-1072 Guide RNA for Insulin

24 samples per tube

Kit contents	Part number	Quantity	Storage temperature
Guide RNA for Insulin	REA1054	1	-80°C

Required materials and equipment provided by the user

Kit contents	Part number	Quantity
Plastics	0.5 mL, 1.5 mL DNA LoBind Tubes	Eppendorf 022431021
	0.2 mL thin wall PCR tubes	General Lab Supplier
	Low-Retention Filtered Sterile Tips (10 µl, 20 µl, 200 µl and 1000 µl)	General Lab Supplier
Reagents	cDNA-amplified and adapter-ligated product from Chromium Next GEM Single Cell 3' Reagent Kits protocol v3.1	10X Genomics
	AMPure XP Beads	Beckman Coulter A63881
	Absolute Ethanol, 200 Proof	General Lab Supplier
	Qubit dsDNA HS Assay Kit	ThermoFisher Scientific Q32854
Equipment	Single Channel Pipettes (10 µl, 20 µl, 200 µl, and 1000 µl)	General Lab Supplier
	Multichannel Pipettes (10 µl, 20 µl, and 200 µl)	General Lab Supplier
	Vortex Mixer	General Lab Supplier
	Microcentrifuge	General Lab Supplier
	PCR Magnetic Rack or Stand for use with tubes	General Lab Supplier
	Ice Bucket	General Lab Supplier
	PCR Thermal Cycler	General Lab Supplier
	Qubit Fluorometer	ThermoFisher Scientific Q33238
	Automated electrophoresis such as TapeStation	General Lab Supplier
DNA analysis instrument, such as the Agilent 2100 Bioanalyzer® System	General Lab Supplier	

Best Practices

General:

- Periodically, optimizations and revisions are made to the kit components and user manual. Therefore, it is important to follow the protocol included with or appropriate for the kit in question. This can be done by comparing the name and version number of the Jumpcode product to the name and version number of the protocol. If you need further assistance in this regard, contact support@jumpcodegenomics.com.
- This protocol describes the reagents, best practices, workflow and method details for DepleteX™ depletion as it applies to the 10x Genomics Chromium Next GEM Single Cell 3' protocol v3.1. Please refer to the Chromium Next GEM Single Cell 3' protocol for warnings and precautions related to Chromium library preparation reagents.
- Use properly calibrated pipettes as library preparation is sensitive to pipetting error.
- Ensure that all pipette tips, microcentrifuge tubes, and other consumables are DNase- and RNase-free.
- Follow good laboratory practices throughout the protocol, including RNA-safe procedures when handling RNA.

Reagent Handling

- Do not remove Cas9 and RNase Inhibitor from -20°C until before use. Return to -20°C immediately after use.
- Store the Guide RNA at -80°C. Do not remove the reagent from -80°C until time of use. Return it to -80°C immediately after use.
- We recommend a maximum of 3 freeze-thaw cycles for the Guide RNA. The Guide RNA tube contains material for 24 samples. It is strongly recommended that multiple smaller aliquots of the Guide RNA be prepared when the reagent is first thawed in order to reduce the number of freeze-thaw cycles affecting the Guide RNA.
- Do not freeze AMPure® XP beads.
- Allow AMPure XP beads to come to room temperature for 30 minutes before use.
- Vortex AMPure XP beads immediately before use. Ensure that they are in a uniform suspension before use.
- Use magnetic stands appropriate for PCR tubes

Revision log

Version	Date	Description
V1.0	September 2023	Launch Release

Protocol

Before starting DepleteX™ Single Cell Boost Kit protocol

1. Follow the Chromium Next GEM Single Cell 3' protocol v3.1 up to and including Step 3.4.
2. At the end of **Step 3.4: Post Ligation Cleanup – SPRIselect**, elute the library in 16 µL of Nuclease-Free Water or Buffer EB.
3. Transfer 15 µL of the sample to a new 0.2 mL or 0.5 mL nuclease-free microcentrifuge tube.
4. Proceed to the **Step A: Ribonucleoprotein complex formation**.

Step A: Ribonucleoprotein complex formation

 **Hands-on time: 5 min | Total time: 15 min**

Reagents preparation

Item	Storage	Handling
10X Cas9 Buffer	-20°C	If frozen, thaw on ice. Finger-tap mix and spin down. Keep on ice. Return to freezer immediately after use.
RNase Inhibitor		
Cas9		
Guide RNAs (2 tubes)	-80°C	

The first step in the protocol is split into sub-sections (Part A and Part B) based on variations in use of different products.

If using only one set of guide RNA, then follow sub-section instructions for **Part A** below:

Each of the following is considered as one set of guide RNA when used by itself.

- KIT1018 Single Cell RNA Boost Kit - Boost guides only (no non-variable genes)
- KIT1024 RNA Depletion Panel (Globin)
- KIT1025 RNA Depletion Panel (Insulin)

If using two sets of guide RNA together, then follow sub-section instructions for **Part B** below:

- KIT1018 DepleteX™ Single Cell RNA Boost Kit - Boost AND non-variable gene depletion
- KIT1018 DepleteX™ Single Cell RNA Boost Kit in combination with either:
 - KIT1024 RNA Depletion Panel (Globin)
 - KIT1025 RNA Depletion Panel (Insulin)

Part A

1. If using one set of Guide RNA by itself, then combine the following reagents in the order listed below in a 0.2 mL or 0.5 mL nuclease-free microcentrifuge tube.

"RNP Complex Formation" Reaction Mix for using one set of Guide RNA		
	Single Cell RNA Boost (KIT1018) - Boost only, no NVG	Globin (KIT1024) or Insulin (KIT1025)
10X Cas9 Buffer	1.0 μ L	1.0 μ L
RNase Inhibitor	1.0 μ L	1.0 μ L
Guide RNA	3.5 μ L	1.6 μ L
Cas9	2.3 μ L	0.9 μ L
Nuclease-free water	1 μ L	4.3 μ L
Total Volume	8.8 μL	8.8 μL


Part B

1. If using a **combination of Guide RNA sets**, then combine the following reagents in the order listed below in a 0.2 mL or 0.5 mL nuclease-free microcentrifuge tube.

"RNP Complex Formation" Reaction Mix for using a combination of Guide RNA sets		
	DepleteX Single Cell RNA Boost (KIT1018) - Boost + NVG	Single Cell RNA Boost combined with either Globin or Insulin
10X Cas9 Buffer	1.0 μ L	1.0 μ L
RNase Inhibitor	1.0 μ L	1.0 μ L
Guide RNA for Single Cell	3.5 μ L Boost	3.9 μ L
Guide RNA set 2	1.0 μ L NVG	1.6 μ L
Cas9	2.3 μ L	3.2 μ L
Total Volume	8.8 μL	10.7 μL

2. Mix the contents gently by pipetting up and down. Centrifuge briefly to collect the contents at the bottom of the tube.
3. Leave the tube on the laboratory bench for 10 minutes at room temperature. This is the ribonucleoprotein complex (RNP).
4. Proceed immediately to **Step B: CRISPR digestion**.

Step B: CRISPR digestion

 **Hands-on time: 5 min | Total time: 65 min**

Reagents preparation

Item	Storage	Handling
10X Cas9 Buffer	-20°C	If frozen, thaw on ice. Finger-tap mix and spin down. Keep on ice. Return to freezer immediately after use.
Ribonucleoprotein complex (RNP) from Step A		
Adapter-ligated cDNA library from Step 3.4 of the Chromium Next GEM Single Cell 3' Reagent Kits protocol v3.1		
AMPure XP beads	4°C	Thaw the AMPure XP beads on the laboratory bench to bring them to room temperature. A minimum of 30 minutes at room temperature is recommended before use.

The first step in the protocol is split into sub-sections (Part A and Part B) based on variations in use of different products. See the beginning of the protocol Step A for definitions of Part A and Part B.

Part A

1. If using **one set of Guide RNA by itself**, then combine the following reagents in a 0.5 mL nuclease-free microcentrifuge tube:

"CRISPR digestion" Reaction Mix for using one Guide RNA set	Volume
10X Cas9 Buffer	1.5 µL
Adapter-ligated cDNA library from Step 3.4 of the Chromium Next GEM Single Cell 3' protocol	15 µL
Ribonucleoprotein complex (RNP) from Step A	8.8 µL
Total Volume	~25 µL

Part B

1. If using a **combination of Guide RNA sets**, then follow the appropriate table below to combine reagents in the order listed below in a 0.5 mL nuclease-free microcentrifuge tube.

"CRISPR digestion" Reaction Mix for using a combination of Guide RNA sets			
	DepleteX Single Cell RNA Boost with NVG	DepleteX Single Cell RNA Boost in combination with Globin or Insulin	
Nuclease-free water	n/a		2.3 μ L
10X Cas9 Buffer	1.5 μ L		2.0 μ L
Adapter-ligated cDNA library from Step 3.4 of the Chromium Next GEM Single Cell 3' protocol	15 μ L		15 μ L
Ribonucleoprotein complex (RNP) from Step A	8.8 μ L		10.7 μ L
Total Volume	~25 μL		~30 μL

2. Mix the CRISPR digestion reaction gently by pipetting up and down. Centrifuge briefly to collect the contents at the bottom of the tube. Incubate the tube at 42°C for 60 minutes.

Note: During this incubation period, remove the AMPure XP beads from the refrigerator and place them on the laboratory bench to bring them to room temperature. A minimum of 30 minutes at room temperature is recommended before use.

3. After the incubation is complete, transfer the tube to ice for ~2 minutes.
4. Briefly centrifuge the tube to collect the contents at the bottom of the tube. Transfer the tube to the laboratory bench. Immediately proceed to **Step C: Size Selection**.

Step C: Size Selection

 **Hands-on time: 15 min | Total time: 45 min**

Reagents preparation

Item	Storage	Handling
Nuclease-Free Water	-20°C	If frozen, thaw on ice. Finger-tap mix and spin down.
CRISPR-digested product from Step B		Keep on ice. Return to freezer immediately after use.
AMPure XP beads	4°C	Thaw the AMPure XP beads on the laboratory bench to bring them to room temperature. A minimum of 30 minutes at room temperature is recommended before use.
80% Ethanol	Room temperature	Freshly prepared

The first step in the protocol is split into sub-sections (Part A and Part B) based on variations in use of different products. See the beginning of the protocol Step A for definitions of Part A and Part B.

- Part A:** If using each **Guide RNA set by itself**, then add 25.3 μL of Nuclease-Free Water to the CRISPR digestion reaction.

Part B: If using a **combination of Guide RNA sets**, then add 20 μL of Nuclease-Free Water to the CRISPR digestion reaction.
- Mix gently by pipetting up and down several times. Place the tube on the laboratory bench at room temperature.
- Add 30 μL of well-resuspended, room-temperature AMPure XP beads, pipette up and down several times to mix and incubate the tube at room temperature for 10 minutes. Mix the sample halfway through the incubation by pipetting up and down several times.
- Place the tube on the magnetic stand. Allow the solution to clear (3–5 minutes). Remove and discard the supernatant without disturbing the beads.
- Add 200 μL of freshly prepared 80% ethanol to the tube.
- After 30 seconds, remove and discard the ethanol. Leave the tube on the magnetic stand during this step.
- Repeat the wash step (previous two steps) with 200 μL of 80% ethanol. Remove as much ethanol as possible after the second wash.
- Open the cap of the tube while it is on the magnetic stand and allow the beads to air dry for 5 minutes.
- Add 31 μL of Nuclease-Free Water to the beads.
- Remove the tube from the magnetic stand. Mix the contents by pipetting up and down several times to fully resuspend the beads in the liquid.
- Incubate the tube at room temperature for 10 minutes. Mix the sample halfway through the incubation (5 minutes) by pipetting up and down several times.
- Place the tube on the magnetic stand. Allow the solution to clear.

- Transfer 30 μL of the supernatant to a new microcentrifuge tube. Place the tube on ice. This tube contains the DepleteX™-depleted Chromium library.
- Proceed to the beginning of **Step 3.5: Sample Index PCR** of the Chromium Next GEM Single Cell 3' protocol v3.1. **Before proceeding, please read the following notes regarding Steps 3.5 and 3.6.**

Step 3.5: Sample Index PCR: Refer to recommendations in the Chromium Next GEM Single Cell 3' protocol to determine the ideal number of PCR cycles. No adjustment in PCR cycles is necessary to compensate for DepleteX™ depletion.

Step 3.6: Post Sample Index PCR Double Sided Size Selection – SPRIselect: Elute the library in 20 μL of Buffer EB, instead of 35.5 μL of Buffer EB, to increase the concentration of the DNA library.

Library validation

It is recommended that the user assess library yield using a dsDNA-specific fluorescence-based method (such as a Qubit fluorometer) and library fragment profile on an Agilent Bioanalyzer 2100 or equivalent instrument before sequencing. Depleted libraries have a similar fragment profile to those of standard Chromium single cell libraries with a typical shift of 20-50 bp to the left of the electropherogram (i.e., the average fragment size is 20-50 bp smaller than that of a standard Chromium library).

It is also recommended that qPCR quantification be performed to ensure optimal cluster density on an Illumina sequencing instrument.

Once the library has been quantitated, it is ready for cluster generation on an Illumina instrument. Please follow standard Illumina protocols for the loading of the library and for cluster generation on the instrument. If the library needs to be stored before sequencing, please store it at -20°C .

Bioinformatics guidelines

10x Genomics Chromium single cell data generated with the DepleteX Single Cell RNA Boost Kit can be analyzed using the Cell Ranger pipeline and other commonly used single cell data analysis tools.

How to compare DepleteX-depleted data to the control data

The standard data analysis workflow must be modified to observe the impact of DepleteX™ depletion on metrics that are typically used to assess single cell library quality, such as numbers of genes per cell and numbers of UMIs per cell. Additionally, guidelines are provided to remove dead cells from downstream analysis. UMAP visualization is also available in the GitHub site.

Use Cell Ranger pipeline with the standard transcript

Follow Jumpcode GitHub instructions in python or R to remove targets to see genes/cell and UMI/cell benefits

Perform subsequent analysis (coverage benefits, ribo/mito/NVG depletion plots, UMAP plots) with the full data set

Warning: If data from DepleteX-depleted libraries is aligned to the comprehensive standard human Cell Ranger Index to perform a comparative analysis between standard Chromium data and equivalent DepleteX-derived data, no significant improvement in the number of UMIs and genes would be observed with the latter. This is the case because the genes targeted for depletion typically constitute ~35% of all Chromium sequencing reads and a significant proportion of unique reads in the sequencing data. When molecules derived from these genes are removed from the library, so are a significant proportion of UMIs in the library.

Access multiple options to support your analysis needs on our Quick Start page:

<https://www.jumpcodegenomics.com/resources/quick-start/crisprclean-single-cell>

- Video walkthrough with our public dataset files.
- Down sampling guidelines for Cell Ranger
- Step-by-step guide in Python or R on how to analyze your data and see the benefits of depletion
- Generate graphs such as the following:
 - Ribo, Mito, and total depletion including NVG content
 - Genes detected and UMI benefit
 - UMAP plots

This product is intended for research purposes only.

This product is not intended to be used for therapeutic or diagnostic purposes in humans or animals.

This product is covered by one or more patents, trademarks and/or copyrights owned or controlled by Jumpcode Genomics, Inc. For more information about commercial rights, please email us at support@jumpcodegenomics.com. While Jumpcode Genomics develops and validates its products for various applications, the use of this product may require the buyer to obtain additional third party intellectual property rights for certain applications.

AMPURE® is a registered trademark of Beckman Coulter, Inc.

BIOANALYZER® is a registered trademark of Agilent Technologies, Inc.

ILLUMINA® is a registered trademark of Illumina, Inc.

DepleteX™ is a trademark of Jumpcode Genomics, Inc.

10X Genomics®, Cell Ranger®, Next GEM®, and Chromium® are registered trademarks of 10X Genomics, Inc.

Qubit® is a registered trademark of Thermo Fisher Scientific, Inc.